

~~TOC~~

## Pile-Up Effect



The Pile-Up effect describes the effects of photons lost at high photon count rates due to the dead time of the TCSPC devices.

Most single photon counting detectors and TCSPC electronics have a **dead time**. After registering one photon the device needs some time to be ready for the next photon to detect. The dead time for TCSPC devices used for the measurement of fluorescence decays are in the range of 25 to 90 ns.

Let's take as an example a laser repetition rate of 40 MHz. The fluorescence molecules receive every 25 ns a laser pulse. Considering the dead time we can detect only one photon after one laser shot. "Pile up" describes the effects of photons lost at high photon count rates due to the dead time of the TCSPC devices. There are two effects:

- The average measured lifetime becomes shorter
- A mono-exponential decay becomes bi-exponential with an additional shorter component.

Since the pile-up effect distorts the histogram, it must be avoided. To do so, the measured count rate should not exceed a certain limit in respect to the laser repetition rate.

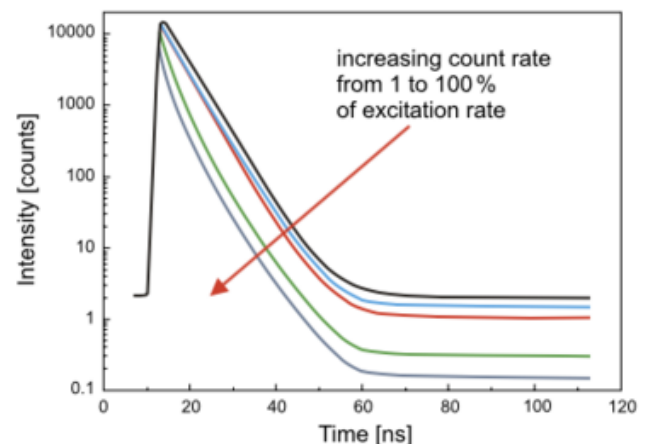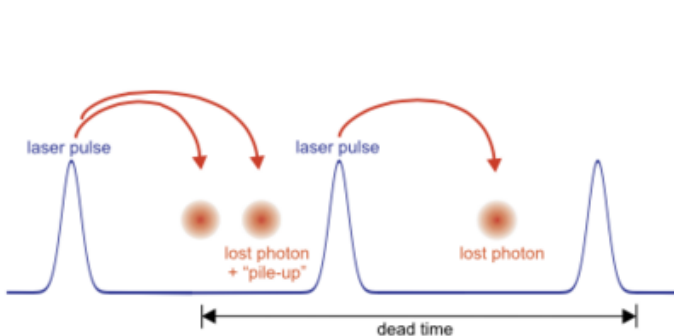

## Practical considerations


For FLIM measurements, the count rate must be no more than 1 - 10% of the excitation rate. This allows the single photon statistics: The probability of detecting more than one fluorescence photon per excitation pulse is negligible.

- [How to Avoid the Pile-up Effect in FLIM Measurements](#)

## Example of pile up effect

We have performed an example measurement with a fluorophore having a lifetime of 4 ns. At 20 MHz laser repetition rate and 2 MHz detection count rate (10%) and a total count number of 10.000 photons in the TCSPC histogram the following results are obtained:

- the average lifetime decrease is 3%
- the  $\chi^2$  value increases from 1.08 to 1.10
- an additional short component lifetime could still not be detected

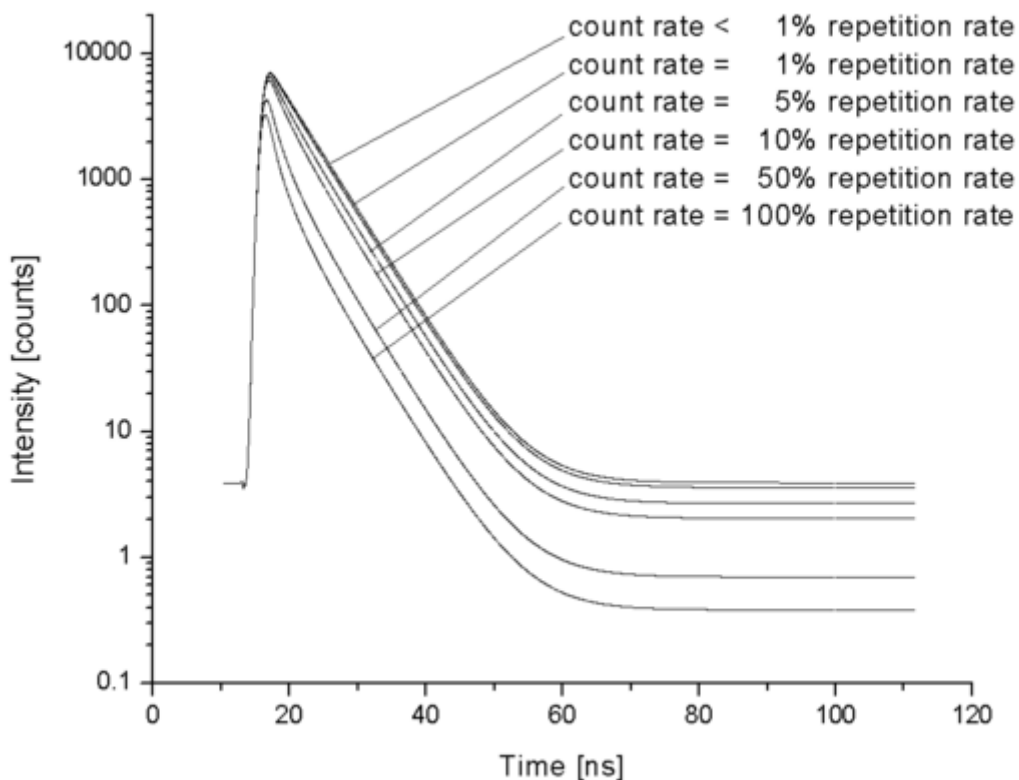

Copyright of this document belongs to PicoQuant GmbH. No parts of it may be reproduced, translated or transferred to third parties without written permission of PicoQuant GmbH. All information given here is reliable to our best knowledge. However, no responsibility is assumed for possible inaccuracies or omissions. Specifications and external appearances are subject to change without notice.



PicoQuant GmbH  
Rudower Chaussee 29 (IGZ)  
12489 Berlin  
Germany

P +49-(0)30-1208820-89  
F +49-(0)30-1208820-90  
info@picoquant.com  
www.picoquant.com